


Standard Controller (Rectangle Controller)

To View Min/Max 'Air' Temperature



STEP 1
RUNNING LOAD TEMPERATURE
The number shown on a **resting display** is the **running load temperature**



STEP 2
Press and quickly release button 
- "thi" will appear on the display



STEP 3
Press and hold button  again to display the MAXIMUM temperature value

Lec Medical Temperature Monitoring Record

Time	Temp	Comments	Site	RL2	Reset	Comments	Site
10:00	44.1						
10:05	44.1						
10:10	44.1						
10:15	44.1						
10:20	44.1						
10:25	44.1						
10:30	44.1						
10:35	44.1						
10:40	44.1						
10:45	44.1						
10:50	44.1						
10:55	44.1						
11:00	44.1						

STEP 4
Record the temperature
You can download min/max tables from:
www.lec-medical.co.uk/support/min-max-tables



STEP 5
Release button  - "tlo" will appear on the display



STEP 6
Press and hold button  again to display the MINIMUM temperature value



STEP 7
T1 will appear, ignore as this function is not used

Lec Medical Temperature Monitoring Record

Time	Temp	Comments	Site	RL2	Reset	Comments	Site
10:00	44.1						
10:05	44.1						
10:10	44.1						
10:15	44.1						
10:20	44.1						
10:25	44.1						
10:30	44.1						
10:35	44.1						
10:40	44.1						
10:45	44.1						
10:50	44.1						
10:55	44.1						
11:00	44.1						

STEP 8
Record the temperature
You can download min/max tables from:
www.lec-medical.co.uk/support/min-max-tables

To Reset Min/Max 'Air' Temperature



STEP 1

Press button  - "thi" will appear on the display





STEP 2

Press and hold the  button again to display the MAXIMUM temperature value




STEP 3

Keeping  held, press button  at the same time. The maximum has now been reset




STEP 4

Release both buttons - "tlo" appears on the screen. Press and hold button  again to display the MINIMUM temperature value



STEP 5

Keeping  held, press button 



STEP 6

The MINIMUM has been reset and the new values will be stored



STEP 7

T1 will appear, ignore as this function is not used



STEP 8

After 5 seconds the display will return to normal, or press button 